




QUV Spectral Power Distribution for QUV with UVA-340 Fluorescent Lamps

SPD for QUV[®] with UVA-340 Fluorescent Lamps

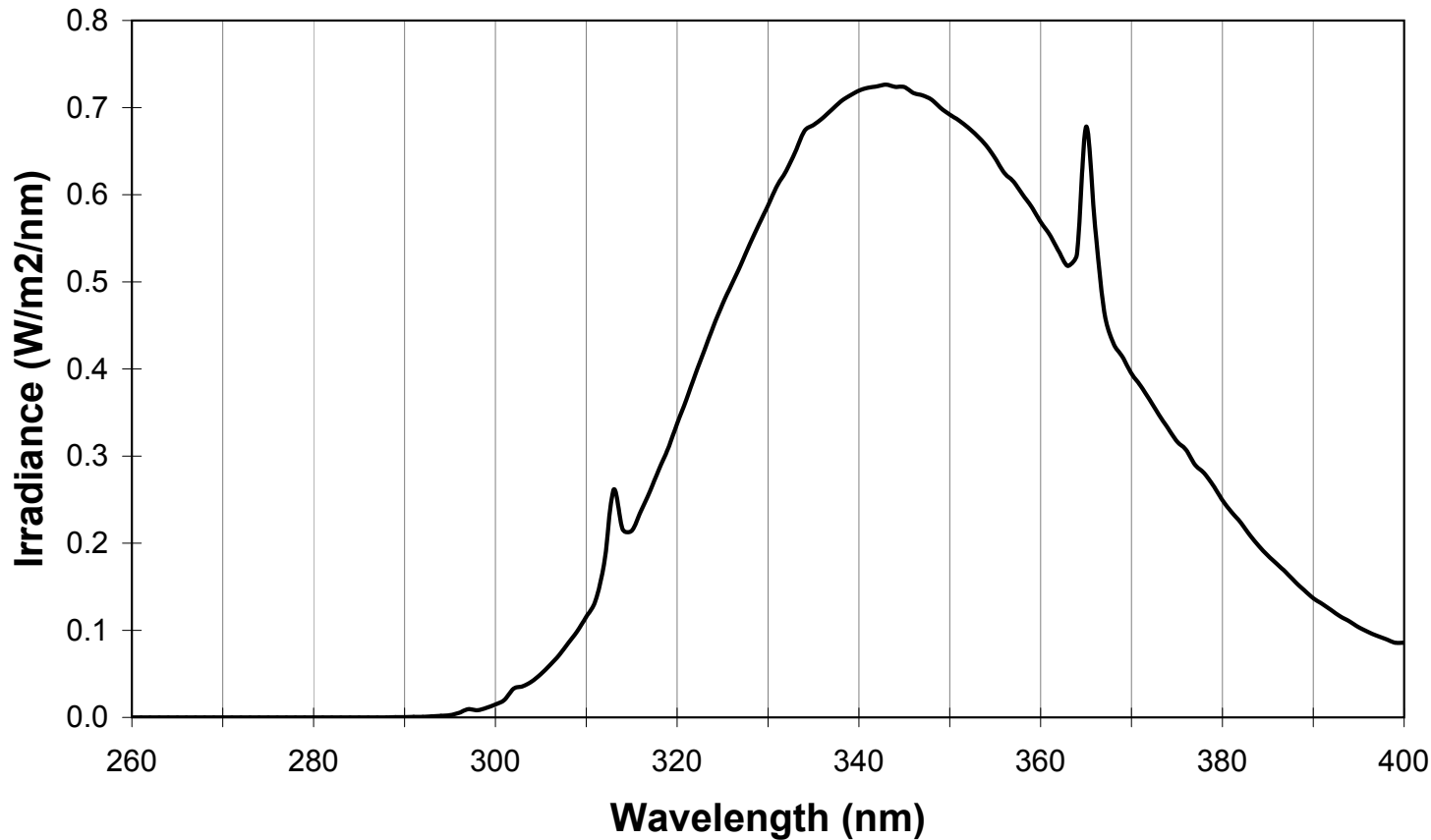

SPD for QUV with UVA-340 Fluorescent Lamps

NOTE: Irradiance values may differ from those listed below due to variations in lamps, lamp age, chamber or room temperature, and measurement position. Different measurement instrumentation and calibration techniques can also cause measured irradiance values to differ from those listed below.

Wave-length (nm)	Irradiance (W/m2/nm)	Wave-length (nm)	Irradiance (W/m2/nm)	Wave-length (nm)	Irradiance (W/m2/nm)	Wave-length (nm)	Irradiance (W/m2/nm)
260	0.0000	295	0.0026	330	0.5877	365	0.6780
261	0.0000	296	0.0052	331	0.6102	366	0.5628
262	0.0000	297	0.0095	332	0.6274	367	0.4653
263	0.0000	298	0.0083	333	0.6491	368	0.4299
264	0.0000	299	0.0110	334	0.6731	369	0.4139
265	0.0000	300	0.0149	335	0.6800	370	0.3947
266	0.0000	301	0.0201	336	0.6879	371	0.3807
267	0.0000	302	0.0330	337	0.6974	372	0.3647
268	0.0000	303	0.0357	338	0.7070	373	0.3476
269	0.0000	304	0.0412	339	0.7137	374	0.3323
270	0.0000	305	0.0496	340	0.7195	375	0.3169
271	0.0000	306	0.0600	341	0.7227	376	0.3073
272	0.0000	307	0.0714	342	0.7244	377	0.2899
273	0.0000	308	0.0850	343	0.7263	378	0.2803
274	0.0000	309	0.0989	344	0.7238	379	0.2662
275	0.0000	310	0.1156	345	0.7236	380	0.2494
276	0.0000	311	0.1339	346	0.7168	381	0.2359
277	0.0000	312	0.1789	347	0.7140	382	0.2238
278	0.0000	313	0.2613	348	0.7091	383	0.2096
279	0.0000	314	0.2160	349	0.6996	384	0.1969
280	0.0000	315	0.2142	350	0.6919	385	0.1859
281	0.0000	316	0.2365	351	0.6854	386	0.1761
282	0.0000	317	0.2594	352	0.6772	387	0.1665
283	0.0000	318	0.2846	353	0.6678	388	0.1558
284	0.0000	319	0.3086	354	0.6566	389	0.1462
285	0.0000	320	0.3371	355	0.6417	390	0.1369
286	0.0000	321	0.3644	356	0.6250	391	0.1303
287	0.0000	322	0.3935	357	0.6151	392	0.1230
288	0.0000	323	0.4215	358	0.6005	393	0.1158
289	0.0003	324	0.4490	359	0.5860	394	0.1102
290	0.0003	325	0.4745	360	0.5688	395	0.1036
291	0.0006	326	0.4969	361	0.5539	396	0.0982
292	0.0006	327	0.5195	362	0.5351	397	0.0937
293	0.0012	328	0.5433	363	0.5183	398	0.0899
294	0.0018	329	0.5659	364	0.5317	399	0.0858
						400	0.0858

Q-Lab Corporation

www.q-lab.com



Q-Lab Headquarters
Westlake, OH USA
Tel: +1-440-835-8700
info@q-lab.com

Q-Lab Florida
Homestead, FL USA
Tel: +1-305-245-5600
q-lab@q-lab.com

Q-Lab Europe, Ltd.
Bolton, England
Tel: +44-1204-861616
info.eu@q-lab.com

Q-Lab Arizona
Buckeye, AZ USA
Tel: +1-623-386-5140
q-lab@q-lab.com

Q-Lab Deutschland, GmbH
Saarbrücken, Germany
Tel: +49-681-857470
vertrieb@q-lab.com

Q-Lab China 中国代表处
Shanghai, China 中国上海
电话: +86-21-5879-7970
info.cn@q-lab.com